


Stirring the Liquids

Before connecting the power supply, make sure that the local mains power voltage matches the stirrer requirements.

1. Place a stir bar into the container.
2. Rotate the control knob clockwise to turn on the stirrer and adjust the stirring speed.
3. If the stir bar is not running smoothly, move container slightly, the stir bar should be in the center of the container.

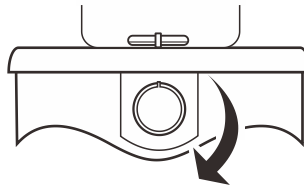

Specification

Maximum Stirring Volume	2000 ml (H ₂ O)
Stirring Speed	0 to 1250 rpm
Top Plate Diameter	145 mm (5.7 in.)
Material	PMMA
Stir Bar Size	30 (L) × 7 (Ø) mm (1.18 × 0.27 in.)
Power Requirements	AC 220V/50Hz
Dimensions	185 (Ø) × 75 (H) mm (7.28 × 2.95 in.)
Weight	0.6 kg (1.3 lb)

Disposal

This product is required to comply with the European Union's Waste Electrical and Electronic Equipment (WEEE) Directive 2002/96/EC and may not be disposed of in domestic waste. Please dispose of product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.



Warranty

The warranty period for stirrer is one year from the date of shipment. Above warranty does not cover the stir bar. Out of warranty products will be repaired on a charged basis. The warranty on your stirrer shall not apply to defects resulting from:

- Improper or inadequate maintenance by customer
- Unauthorized modification or misuse
- Operation outside of the environment specifications of the products

For more information, please contact the supplier.